BFI pharmaceutical



I Application

The range of BFI series presents vertical direct motor-driven agitators.

These agitators are used in processes of mixing, dissolution, dispersion and maintenance whenever powerful and high-speed agitation is required in the food processing, cosmetics, pharmaceutical or chemical industries.

The agitators are normally fitted to small or medium volume tanks for work with low-viscosity products.

Bearing support of the BFI series agitators is manufactured in stainless steel and it is indicated for the pharmaceutical applications.

I Operating principle

The agitators are normally installed inside the tank either centrally with current deflectors, or offset so that the offsetting itself prevents the rotation of the stirred product.

The rotation of the propeller makes the fluid flow to the bottom and then up to the surface of the product along the walls of the tank. The effect is promoted if the bottom of the tank is curved.

The saw tooth propeller promotes the dispersion of the product.



I Design and features

Vertical agitator.

Double lip seal.

Double ball bearing support.

The propeller fixed to the shaft by means of a threaded connection and a gasket; the shaft is made of one piece.

IEC motors, 1500 rpm, IP 55, F-class insulation.

Maximum power 5,5 kW.

Marine propeller (Type 10) or saw tooth propeller (Type 4). Indicated for work under pressure or in vacuum (-1 to 4 bar).



Parts in contact with the product

Retainer

Bearing support

Surface finish

AISI 316

GYLON according to FDA 177.1550

AISI 316

Electropolished





BFI pharmaceutical

I Options

Double/flushed mechanical seal.


Shaft and propeller lined with plastic material for use with corrosive products. Welded propeller with Ra $\leq 0.4~\mu m$ surface finish.

Motor shroud.

Motors with other protections.

Motors: 950 rpm or 750 rpm.

Geared motors or variable speed drives.

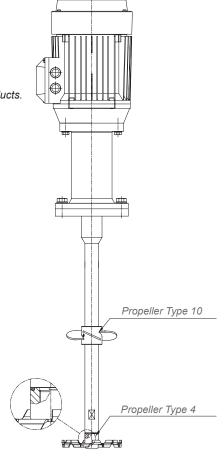

I Technical specifications and dimensions

			ize							Туре		
	e									10	4	
Agitator type	Motor power	Speed Bearing support size		Dimensions					Agitator shaft		Marine propeller	Cowler
	FLAME	[rpm]	Be	С	Flange				a al		øΗ	~ 11
	[kW]				ø A	ø B	DN	ø G(n)	ø d	L _{max} .	ØH	øΗ
BFI 1.10-4011-1-160	1.1	1450	90	515				14(4)	20	1400	160	
BFI 1.10-4015-1-175	1.5										175	
BFI 1.10-4022-1-200	2.2		100	543	210	180	100				200	
BFI 1.10-4030-1-200	3										200	
BFI 1.10-4040-1-225	4		112	546							225	
BFI 1.10-4055-1-225	5.5										225	
BFI 1.10-6007-1-175	0.75	950	90	515							175	
BFI 1.10-6011-1-200	1.1										200	
BFI 1.10-6015-1-225	1.5		100	543							225	
BFI 1.10-6030-1-250	3		112	546							250	
BFI 1.4-4022-1-150	2.2	1450	100	543					45	1500		150
BFI 1.4-4040-1-200	4	1430	112	546								200







